

Who's Guarding You & Your Animals?

Peroxigard® Concentrate One-Step Disinfectant Cleaner for use in Life Sciences

One technology for all your facility's cleaning and disinfection needs.

Efficacy

Effective against over 75+ key pathogens, including those on bioexclusion lists. Achieves disinfection up to 2x faster than other concentrate products.

Safety

Designed with animal safety in mind. Fragrance free, non-irritating to eyes and skin, non-corrosive and superior compatibility at the in-use dilution. Leaves no active residue on surfaces or equipment.

Simplicity

Easy to dilute concentrate, with instant activation. Diluted solution lasts up to 90 days, significantly longer than unstable chlorine dioxides. Saving you and your staff time.



Disinfection Compliance for Better Science

Peroxigard Concentrate utilizes Accelerated Hydrogen Peroxide®, a patented and synergistic blend of commonly used ingredients, that when combined with low levels of hydrogen peroxide create an ideal one-step cleaner and disinfectant. Trusted EPA Registered disinfectant with proven efficacy claims and excellent safety profile.

Peroxigard Concentrate is a simple and effective way to clean and disinfect floors, ceilings and walls within your facility. Peroxigard Concentrate is proven to remain effective in the presence of a 5% soil load, providing disinfection confidence. Ideal for use with a variety of equipment, including mixing stations, application equipment/foamers and more.

Effective against key pathogens in as little as 5 minutes* including:

Bacteria

Klebsiella pneumoniae; *Methicillin-resistant Staphylococcus aureus* (MRSA); *Salmonella enterica*; *Mycoplasma orale*

Viruses

Murine Parvovirus; *HIV Type 1*; *Rotavirus*; *Poliovirus Type 1*; *Canine Parvovirus*

Fungi

Microsporium canis; *Trichophyton interdigitale*



DILUTION MADE EASY

Ratio	Per Gallon of Water	Per 4 Gallons of Water
1:16	8.0 fl.oz.	32.0 fl.oz.
1:64	2.0 fl.oz.	8.0 fl.oz.

FOR USE IN:

- Vivariums
- Breeding facilities
- Lab Animal Research facilities
- Life Science laboratories
- Animal Examination rooms
- Non-human primate facilities

FOR USE ON:

- Floors, ceilings, and walls
- Lab benches and exam tables
- Cages, crates, and carts
- Animal housing areas

SHELF LIFE

- 3 years (Opened/Unopened)
Remains effective for up to 90 days once diluted

AVAILABLE FORMATS

- 1 gal, 5 gal, 55 gal
(EPA Reg No. 74559-4)

To learn more about Accelerated Hydrogen Peroxide or check out our peer-reviewed studies, please visit: Peroxigard.com



Peroxigard® is a member of the Virox family of brands.

DISCLAIMER: You must read and follow the instructions on the label for our product and use the product for the particular applications specified on the product label. To the full extent permitted under applicable law, Virox Technologies Inc. shall not be liable for any damages whatsoever, arising out of or otherwise in connection with any failure to follow those instructions or any use of its products which is not specified on the product's label. The instructions on the product label are not a warranty, express or implied, and Virox Technologies, Inc. expressly disclaims any implied warranties, including any warranty of fitness for a particular purpose. Under applicable law, some of the above waivers or disclaimers of warranty, or exclusions of liability for certain types of damages may not apply to you. In such cases, our liability will be limited to the fullest extent permitted by applicable law. Product shots have been digitally altered and are not accurate as shown. Refer to product label or reference sheet for a complete list of disinfection claims and additional features. *See reference sheet for a complete list of efficacy claims, additional features and directions for use.

AHP® and Design and Accelerated Hydrogen Peroxide® and Design are trademarks of Diversey, Inc.

B-091624-8492-08F © 2024 Virox Technologies Inc.